

EzDecap™
Automated Screw Cap Decapper
BRDC-0000

Operation Manual

Ver. 2.0



Copyright © 2024 Blue-Ray Biotech Corp.
All Rights Reserved.

Table of Contents

1. Safety Precautions	1
1.1 Intended Use.....	1
1.2 General Safety.....	1
2. Product Overview	2
3. Installation	2
4. Operation Modes	2
5. Decapping / Capping	3
6. Cycle Time Setting	4
7. Transportation and Storage	5
8. Maintenance	6
8.1 Cleaning and Decontaminating	6
8.2 Gripper Replacement.....	6
9. Disposal	6
Appendix A: Technical Specifications	7
Appendix B: CE Declaration.....	8
Appendix C: Order Information	9

1. Safety Precautions

Before using the **EzDecap™** for the first time, please read this entire Operation Manual carefully. To guarantee problem-free, safe operation of the **EzDecap™**, it is essential to observe the following sections.

1.1 Intended Use

The **EzDecap™** is intended to be used by trained personnel to perform screw cap decapping/capping. In this manual, we assume that the user has knowledge of basic laboratory procedures.

1.2 General Safety



PHYSICAL INJURY HAZARD


Use of the instrument in a manner, or for purposes not specified by Blue-Ray Biotech, may result in personal injury or damage to the instrument.

Using the device in a manner not specified by Blue-Ray Biotech may result in personal injury or damage to the device.

- Do not use the device in a potentially explosive environment or with potentially explosive chemicals.
- Avoid placing the device in direct sunlight.
- Install the device in a location free of excessive dust.
- Install the device in a room with a temperature of 15 - 30 °C, relative humidity of 20 - 80%.
- Choose a flat, stable surface capable of bearing the weight of the device.
- Make sure the power source conforms to the required power supply specifications.
- To avoid electric shock, make sure the device is plugged into a grounded electrical outlet.
- Do not allow water or any foreign objects to enter the various openings of the device.
- The use of excessive force should be avoided.
- Only use the original power adapter supplied by the manufacturer.
- To ensure proper ventilation, make sure the device has at least 30 cm of free space all around.

- Do not place this device where it is difficult to disconnect the power adapter.

2. Product Overview

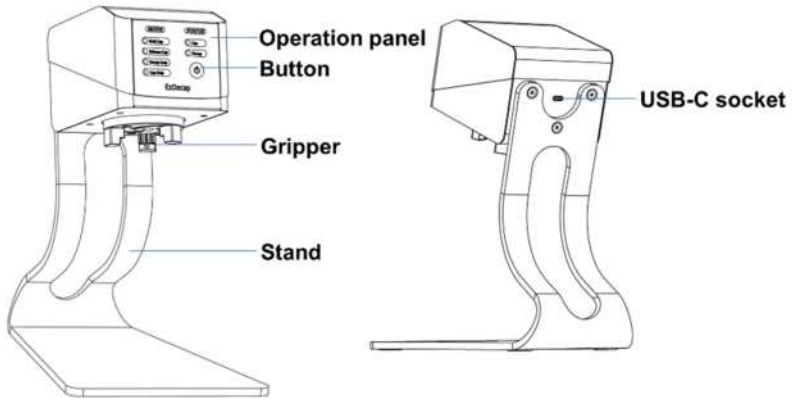




3. Installation

1. Plug the USB-C jack of the power adapter into the USB-C socket on the back of the device.
2. Plug the power adapter into an external power source.

4. Operation Modes

Five operation modes satisfy users with various needs for decapping and capping screw caps.

Indicator	Mode	Action
 Hold Cap	Hold Cap	Decapping with cap held → Capping the held cap
 Release Cap	Release Cap	Decapping with cap released → Capping the cap on the top of the tube
 Decap Only	Decap Only	Decapping and releasing cap immediately

Indicator	Mode	Action
● Cap Only	Cap Only	Capping the cap on the top of the tube
● Release Cap ● Decap Only	Decap and Release	Decapping and releasing cap after 1.5 sec

5. Decapping / Capping

1. Press the power button for 3 seconds to turn on the device.
2. Press the power button to select the mode.
3. If the cap is too tight, it may need to be manually loosened firstly.
4. Hold the tube in hand, and lightly press straight up on the center of the disc to trigger the motor. Once the motor is triggered, you don't need to apply upward force.

Tip I: It could be easier to open the cap by holding the tube and turning it in the opposite direction when the motor is activated.

Tip II: If the cap is not closed correctly, repeat steps 3 and 4.



5. If the cap needs to be extremely tight, it may need to be tightened manually.
6. Press the power button for 3 seconds to turn off the device.

6. Cycle Time Setting

The **EzDecap™** is built with 3 different cycle times for selection to accommodate most tubes. “Shorter time” setting provides express operation, while “Longer time” setting is for the tubes needing longer cycle time. Follow the procedures to set the cycle time.



1. Press the power button for 3 seconds to turn on the device.
2. In “Hold Cap” mode, use a tube to press up on the center of the disc to trigger the motor. The indicators will show: “Hold Cap” in mode and “Cap” in status.



3. Press the power button for 3 seconds to select the different cycle time. The LEDs will blink when setting the cycle time.

Setting	Decapping Cycle Time (sec)	Capping Cycle Time (sec)
Setting 1- Default	1.1	1.0



Setting	Decapping Cycle Time (sec)	Capping Cycle Time (sec)
Setting 2- Shorter time	1.1	0.7
		
Setting 3- Longer time	1.9	1.8
		

- Use a tube again to press up on the center of the disc to trigger the motor and complete the setting.

Note
The cycle time setting will be kept after rebooting the EzDecap™ .

7. Transportation and Storage

This device should be transported and stored in an environment with a temperature between -10 and 60 °C and with a relative humidity of 20 - 80%


8. Maintenance

8.1 Cleaning and Decontaminating

Before cleaning or decontamination, disconnect the power adapter. The decapper can be cleaned or decontaminated with 70% ethanol or a 10% commercial bleach solution.

8.2 Gripper Replacement

The grippers will require occasional replacement, particularly when the grippers can't grab the cap firmly. Use an M2 Allen key to remove the hexagon socket screws on the inside of the grippers. Replace the three grippers with new ones.

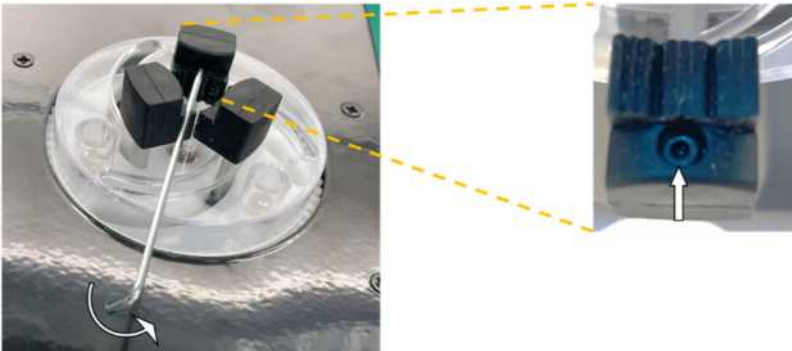

9. Disposal



Do not dispose of this product as unsorted municipal waste.

Follow local municipal waste ordinances for proper disposal provisions to reduce the environmental impact of waste electrical and electronic equipment (WEEE).

European Union Customers: Call your local Blue-Ray Biotech Distributor's Customer Service office for device pick-up and recycling.

Appendix A: Technical Specifications

Dimensions (W x D x H)	170 x 240 x 300 mm
Weight	3.8 kg
Cycle Time Decap/Cap	< 3 sec
Compatibility	11 - 40 mm diameter screw cap tubes
Sensing Method	Pressure sensor
Operation Modes	Five modes
Power Adapter	Input: AC100 - 240 V, 50/60 Hz; Output: DC 24 V, 2.7 A, 65 W
Power Jack	USB-C
Certification	CE

Appendix B: CE Declaration



BLUE-RAY BIOTECH CORP.
5F., No. 2, Aly. 2, Siwei Ln., Zhongzheng Rd.,
Xindian Dist., New Taipei City 231,
Taiwan (R.O.C.)

Declaration of Conformity

Product Name: EzDecap™

Model Names: BRDC-0000

All models comply with the following European standards:

EMC:

EN 55011, BS EN 55011, CISPR 11,

EN IEC 61326-1:2013

EN IEC 61000-3-2:2014

EN 61000-3-3:2013+A1:201

Safety:

EN 61010-1:2010/A1:2019, IEC 61010-1:2010,

IEC 61010-1:2010/AMD1:2016,

BS EN 61010-1:2010+A1:2019

To the best of my knowledge and belief, these units conform to these standards.

Name: **Jimmy Kuo**

Position: Quality Assurance Manager

Issue Date: 2024. 4. 19

Appendix C: Order Information

Order No.	Description
BRDC-0000	EzDecap™ Automated Screw Cap Decapper
BRDC-a001	Gripper set



For more information,
visit www.blue-raybio.com
EzDecap™ Product Webpage

Blue-Ray Biotech Corp.

5F., No. 2, Aly. 2, Siwei Ln., Zhongzheng Rd., Xindian Dist.,
New Taipei City 231, Taiwan (R.O.C.)

P/N: 401-DCBR00-20